



The New Secura®: Easy, Reliable Weighing in Regulated Areas

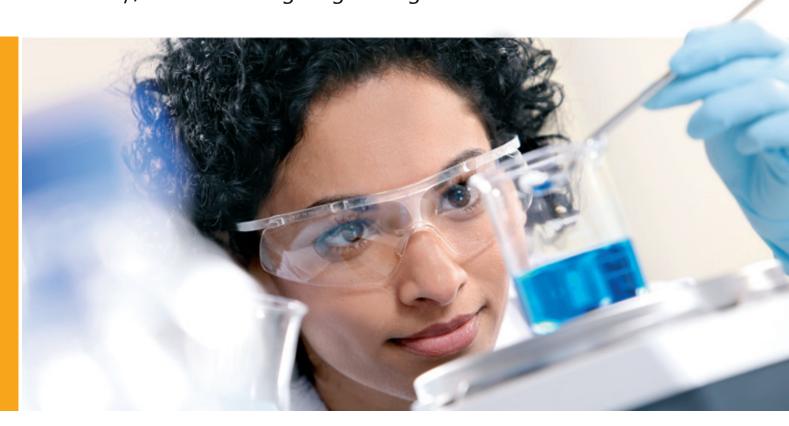







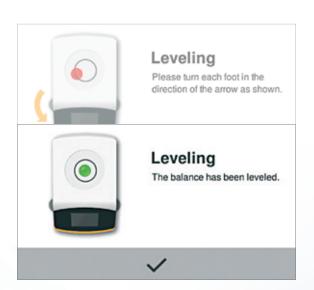

Always Correctly Leveled with Sartorius LevelControl

Secura® has an optoelectronic sensor that continuously checks to see if the balance is perfectly level. If it detects any out-of-tolerances, you will be immediately alerted to this pending risk and see specific guidance prompts to help you level the balance. Output of the weighing data over the interface port will be temporarily interrupted until the balance has been correctly leveled again.

This means you can rest assured knowing that only correct data is transmitted. During leveling, you will not have to change your seated position. More importantly, when working with toxic substances under a laboratory fume hood, you will not have to lean over the balance while leveling it. As a result, Secura rules out the risk of health hazards.

Regulatory Compliance with Expert Support

Do you use your Secura® or other instruments in a regulated area and have to comply with a multitude of requirements, regulations and directives? Relax. Sartorius Lab Products & Services will provide you with the appropriate expert services specifically tailored to ensure compliance with the strict requirements of the pharmaceutical industry – and not just for balances.



Risk Minimization – IQ | OQ Documentation

Do you need qualification documents as proof of suitability to be able to use your Secura® in a regulated area? Sartorius Lab Products & Services will support you by providing trained specialists on site. We

will generate the documents you need according to riskbased test plans and test reports for installation and operational qualification.


Stress-Free USP Minimum Sample Weight Measurement

Compliance with the USP minimum sample weight requirements with absolute process reliability? No problem for the new Secura®.

The display clearly indicates when the minimum sample weight is below the required minimum limit. In addition, the balance will temporarily block data transfer to a printer or a device connected to its interface port. This gives you

peace of mind knowing that inadequate data cannot be inadvertently processed further.

Our specially trained service staff will determine the minimum sample weight for you right at the balance's site of installation, program it into your Secura® and attach a corresponding verification seal to the balance.

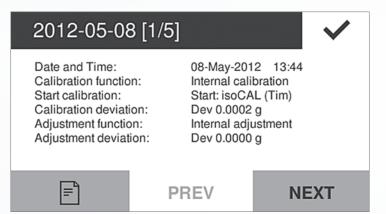

Best Results – Quality Documentation Made Easy

Secura® features a Cal Audit Trail function to document the quality of each weigh-in procedure you perform. Every combined calibration and adjustment procedure and each leveling process, including any non-conformities, are recorded by the balance, can be viewed on the touch screen and, if necessary, can be transferred over the balance's interface to a computer or to a Sartorius laboratory printer.




No Risk Due to Temperature Fluctuations

Even the slightest temperature fluctuations have a major influence on the repeatability and quality of the weighing results.

This is why every Secura® comes equipped with the isoCAL internal calibration and adjustment function that ensures consistently accurate results.

Secura® can be configured to prompt you to launch this procedure or to have isoCAL take care of this as soon as an activation threshold is reached.

The Cal Audit Trail function records each isoCAL procedure performed so it can be traced for your quality assurance.

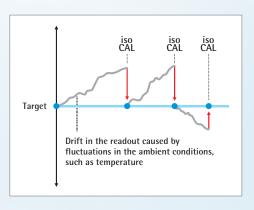

Cleanability to Secure the Best Results

Cleanliness is one of the main prerequisites in the laboratory to prevent contamination and ensure work area safety.

That's why Sartorius placed a premium on designing Secura® to enable easy and thorough cleaning.

Sleek, resistant surfaces plus clear contours and few edges ensure that Secura® will help you stay compliant with all hygiene requirements.





Security by Password Protection

Secura® comes with password protection to restrict adjustment of the balance settings only to authorized persons. If you set a password, all functions that could change the metrological performance specifications are inactivated. This ensures secure process results.

Interactive User Guidance for Fast GLP-Compliant Printouts

If you require traceable documents, Secura® will guide you through the steps for obtaining GLP-compliant printouts. The novel GLP user guidance features will support you, especially for working with traceable sample identification.

Used in conjunction with the new premium YDP30 laboratory printer, which is automatically detected by Secura®, the balance generates printouts of the results on standard printer paper or on continuous labels.







Secura® Technical Specifications

	-1S 1102-1S	2102-1S	3102-1S	5102 - 1S	213-1S	313-1S	513 - 1S	124 - 1S	224-15		Model
Readability mg	7-406 14-557-405	14-557-406	14-557-407	14-557-408	14-557-402	14-557-403	14-557-404	14-557-400	14-557-401).	Fisher Scientific Catalog No
Repeatability mg	1,100	2,100	3,100	5,100	210	310	510	120	220	g	Weighing capacity
Linearity mg 0.2 0.2 2 2 2 2 20 20 20 20 Min. sample weight a cocording to USP, typical Temperature limits for verified models with activated isoCAL Typical stabilization time s 2 2 2 1.5 1.5 1.5 1.5 1.5 1.5 1.5 1.5 1.5 Weighing pan size mm Ø 90 Ø 90 Ø 120 Ø 120 Ø 120 Ø 180 Ø 180 Ø 180 Weighing chamber height mm 209 209 209 209 209 − − − Net weight, approx. kg 5.1 5.1 5.1 5.1 5.1 5.1 5.2 5.2 5.2 5.2 Dimensions, Dimensions, Daw H mm (Inches) Dimensions, Charles and Mini USB - Automatic recognition of Sartorius printer models YDP30 or YDP40 - Direct data transfer to Microsoft Windows programs - Choice of SBI and xBPI data transfer protocols APC functions (Advanced Pharma Compliance) APC functions (Advanced Pharma Compliance) Leveling Display Touch screen with Sartorius graphical user interface optimized for users in pharmaceutical laborat Application programs Weighing, Mixing, Components, Statistics, Density, Percentage, Conversion, Unstable Conditions, Components, Statistics, Density, Percentage, C	10	10	10	10	1	1	1	0.1	0.1	mg	Readability
Min. sample weight according to USP, typical Temperature limits for verified models with activated isoCAL Typical stabilization time s 2 2 2 1.5 1.5 1.5 1.5 1.5 1.5 1.5 1.5 1.5 1.5	10	10	10	10	1	1	1	0.1	0.1	mg	Repeatability
Temperature limits for verified models with activated isoCAL Typical stabilization time s 2 2 2 1.5 1.5 1.5 1.5 1.5 1.5 1.5 1.5 1.5 1.5	20	20	20	20	2	2	2	0.2	0.2	mg	Linearity
for verified models with activated isoCAL Typical stabilization time s 2 2 1.5 1.5 1.5 1.5 1.5 1.5 1.5 1.5 Weighing pan size mm Ø 90 Ø 90 Ø 120 Ø 120 Ø 120 Ø 180 Ø 180 Ø 180 Ø 180 Weighing chamber height mm 209 209 209 209 209 Net weight, approx. kg 5.1 5.1 5.1 5.1 5.1 5.1 5.1 5.1 5.2 5.2 5.2 5.2 (ltb) (11.2) (11.2) (11.2) (11.2) (11.2) (11.2) (11.5) (11.5) (11.5) (11.5) Dimensions, 360×216×320 360×216×95 D×W×H mm (Inches) (14.1×8.5×12.6) (14.1×8.5×3.75) Interface Mini USB - Automatic recognition of Sartorius printer models YDP30 or YDP40 - Direct data transfer to Microsoft Windows® programs - Choice of SBI and xBPI data transfer protocols APC functions (Advanced Pharma Compliance) - Fully automatic temperature- and time-controlled internal calibration and adjustment - isoCAL - Temporary blockage of data transfer to a printer or a computer when uncertain weighing results such as when a result is below the USP minimum sample weight limit, the balance is not level or calibration adjustment needs to be performed - Storage of all data of calibration procedures - Cal Audit Trail Leveling Intelligent, optoelectronic leveling sensor with alarm function and interactive user guidance for reliable leveling Display Touch screen with Sartorius graphical user interface optimized for users in pharmaceutical laborat Application programs Weighing, Mixing, Components, Statistics, Density, Percentage, Conversion, Unstable Conditions, Components, Statistics, Density, P	12	12	12	12	1.5	1.5	1.5	0.12	0.12	g	
Weighing pan size mm Ø 90 Ø 90 Ø 120 Ø 120 Ø 120 Ø 180 Ø 180 Ø 180 Weighing chamber height mm 209 209 209 209 209 – – – Net weight, approx. kg 5.1 5.1 5.1 5.1 5.1 5.2 5.2 5.2 (lb) (11.2) (11.2) (11.2) (11.2) (11.2) (11.2) (11.5) (11.5) (11.5) Dimensions, 360×216×320 360×216×95 D×W×H mm (Inches) (14.1×8.5×12.6) (14.1×8.5×3.75) Interface Mini USB - Automatic recognition of Sartorius printer models YDP30 or YDP40 - Direct data transfer to Microsoft Windows® programs - Choice of SBI and xBPI data transfer protocols APC functions (Advanced Pharma Compliance) — Monitoring of compliance with the USP minimum sample weight limits – SQmin - Password protection of set-up settings Compliance) — Fully automatic temperature— and time—controlled internal calibration and adjustment — isoCAL - Temporary blockage of data transfer to a printer or a computer when uncertain weighing results such as when a result is below the USP minimum sample weight limit, the balance is not level or calibration adjustment needs to be performed - Storage of all data of calibration procedures — Cal Audit Trail Leveling Intelligent, optoelectronic leveling sensor with alarm function and interactive user guidance for reliable leveling Display Touch screen with Sartorius graphical user interface optimized for users in pharmaceutical laborat Application programs Weighing, Mixing, Components, Statistics, Density, Percentage, Conversion, Unstable Conditions, Components, Statistics, Density, Percenta		+10°C +30°C							for verified models with		
Weighing chamber height mm 209 209 209 209 209 Net weight, approx. kg 5.1 5.1 5.1 5.1 5.1 5.1 5.2 5.2 5.2 5.2 (Ib) (11.2) (11.2) (11.2) (11.2) (11.2) (11.5) (11.5) (11.5) (11.5) Dimensions, 360×216×320 360×216×95 D×W×H mm (Inches) (14.1×8.5×12.6) (14.1×8.5×3.75) Interface Mini USB - Automatic recognition of Sartorius printer models YDP30 or YDP40 - Direct data transfer to Microsoft Windows® programs - Choice of SBI and xBPI data transfer protocols APC functions (Advanced Pharma Compliance) - Monitoring of compliance with the USP minimum sample weight limits - SQmin - Password protection of set-up settings Compliance) - Fully automatic temperature- and time-controlled internal calibration and adjustment - isoCAL - Temporary blockage of data transfer to a printer or a computer when uncertain weighing results such as when a result is below the USP minimum sample weight limit, the balance is not level or calibration adjustment needs to be performed - Storage of all data of calibration procedures - Cal Audit Trail Leveling Intelligent, optoelectronic leveling sensor with alarm function and interactive user guidance for reliable leveling Display Touch screen with Sartorius graphical user interface optimized for users in pharmaceutical laborat Application programs Weighing, Mixing, Components, Statistics, Density, Percentage, Conversion, Unstable Conditions, Components, Statistics, De	1.5	1.5	1.5	1.5	1.5	1.5	1.5	2	2	S	Typical stabilization time
Net weight, approx. kg 5.1 5.1 5.1 5.1 5.1 5.1 5.1 5.2 5.2 5.2 5.2 (lb) (11.2) (11.2) (11.2) (11.2) (11.2) (11.5) (11.5) (11.5) (11.5) Dimensions, D×W×H mm (Inches) (14.1×8.5×12.6) 360×216×95 Automatic recognition of Sartorius printer models YDP30 or YDP40 Direct data transfer to Microsoft Windows® programs Choice of SBI and xBPI data transfer protocols APC functions (Advanced Pharma Password protection of set-up settings Compliance) Fully automatic temperature- and time-controlled internal calibration and adjustment – isoCAL Temporary blockage of data transfer to a printer or a computer when uncertain weighing results such as when a result is below the USP minimum sample weight limit, the balance is not level or calibration adjustment needs to be performed Storage of all data of calibration procedures – Cal Audit Trail Leveling Intelligent, optoelectronic leveling sensor with alarm function and interactive user guidance for reliable leveling Display Touch screen with Sartorius graphical user interface optimized for users in pharmaceutical laborat Application programs Weighing, Mixing, Components, Statistics, Density, Percentage, Conversion, Unstable Conditions, Components, Statistics, Density, Percentage, Conversion, Unstable Co	Ø 180	Ø 180	Ø 180	Ø 180	Ø 120	Ø 120	Ø 120	Ø 90	Ø 90	mm	Weighing pan size
(lb) (11.2) (11.2) (11.2) (11.2) (11.2) (11.5) (11.5) (11.5) (11.5) Dimensions, D×W×H mm (Inches) (14.1×8.5×12.6) Mini USB - Automatic recognition of Sartorius printer models YDP30 or YDP40 - Direct data transfer to Microsoft Windows® programs - Choice of SBI and xBPI data transfer protocols APC functions (Advanced Pharma Compliance) - Fully automatic temperature- and time-controlled internal calibration and adjustment – isoCAL - Temporary blockage of data transfer to a printer or a computer when uncertain weighing results such as when a result is below the USP minimum sample weight limit, the balance is not level or calibration adjustment needs to be performed - Storage of all data of calibration procedures – Cal Audit Trail Leveling Display Touch screen with Sartorius graphical user interface optimized for users in pharmaceutical laborat Weighing, Mixing, Components, Statistics, Density, Percentage, Conversion, Unstable Conditions, C	_	-	-	-	209	209	209	209	209	mm	Weighing chamber height
D×W×H mm (Inches) (14.1×8.5×12.6) (14.1×8.5×3.75) Interface Mini USB - Automatic recognition of Sartorius printer models YDP30 or YDP40 - Direct data transfer to Microsoft Windows® programs - Choice of SBI and xBPI data transfer protocols APC functions (Advanced Pharma Compliance) - Monitoring of compliance with the USP minimum sample weight limits – SQmin - Password protection of set-up settings Compliance) - Fully automatic temperature- and time-controlled internal calibration and adjustment – isoCAL - Temporary blockage of data transfer to a printer or a computer when uncertain weighing results such as when a result is below the USP minimum sample weight limit, the balance is not level or calibration adjustment needs to be performed - Storage of all data of calibration procedures – Cal Audit Trail Leveling Intelligent, optoelectronic leveling sensor with alarm function and interactive user guidance for reliable leveling Display Touch screen with Sartorius graphical user interface optimized for users in pharmaceutical laborat Application programs Weighing, Mixing, Components, Statistics, Density, Percentage, Conversion, Unstable Conditions, Componen	5.2 (11.5)									-	Net weight, approx.
- Automatic recognition of Sartorius printer models YDP30 or YDP40 - Direct data transfer to Microsoft Windows® programs - Choice of SBI and xBPI data transfer protocols APC functions - Monitoring of compliance with the USP minimum sample weight limits – SQmin - Password protection of set-up settings Compliance) - Fully automatic temperature- and time-controlled internal calibration and adjustment – isoCAL - Temporary blockage of data transfer to a printer or a computer when uncertain weighing results such as when a result is below the USP minimum sample weight limit, the balance is not level or calibration adjustment needs to be performed - Storage of all data of calibration procedures – Cal Audit Trail Leveling Display Display Touch screen with Sartorius graphical user interface optimized for users in pharmaceutical laborat Weighing, Mixing, Components, Statistics, Density, Percentage, Conversion, Unstable Conditions, Components,										•	
(Advanced Pharma Password protection of set-up settings Fully automatic temperature- and time-controlled internal calibration and adjustment – isoCAL Temporary blockage of data transfer to a printer or a computer when uncertain weighing results such as when a result is below the USP minimum sample weight limit, the balance is not level or calibration adjustment needs to be performed Storage of all data of calibration procedures – Cal Audit Trail Leveling Intelligent, optoelectronic leveling sensor with alarm function and interactive user guidance for reliable leveling Display Touch screen with Sartorius graphical user interface optimized for users in pharmaceutical laborat Application programs Weighing, Mixing, Components, Statistics, Density, Percentage, Conversion, Unstable Conditions, Components, Statistics, Density, Percentage, Conversion, Unstable				DP40	 Automatic recognition of Sartorius printer models YDP30 or Y Direct data transfer to Microsoft Windows® programs 						Interface
user guidance for reliable leveling Display Touch screen with Sartorius graphical user interface optimized for users in pharmaceutical laborat Application programs Weighing, Mixing, Components, Statistics, Density, Percentage, Conversion, Unstable Conditions, C	ts are detected,	g results are d	adjustment – tain weighing	ibration and a	d internal cal or a compute sample weigh	gs me-controlle to a printer of SP minimum performed	set-up settin ature- and ti data transfer below the U needs to be	protection of matic temper blockage of nen a result is a adjustment	PasswordFully autoTemporarysuch as wheelibration		(Advanced Pharma
Application programs Weighing, Mixing, Components, Statistics, Density, Percentage, Conversion, Unstable Conditions, C										Leveling	
	itories	Touch screen with Sartorius graphical user interface optimized for users in pharmaceutical laboratories								Display	
	Checkweighing,	Weighing, Mixing, Components, Statistics, Density, Percentage, Conversion, Unstable Conditions, Checkweighing, Peak Hold, Counting									Application programs
Accessories See Secura® list of accessories		See Secura [®] list of accessories						See Secura®		Accessories	



In the United States:

For customer service, call 1-800-766-7000. To fax an order, use 1-800-926-1166. To order online: www.fishersci.com

In Canada:

For customer service, call 1-800-234-7437. To fax an order, use 1-800-463-2996. To order online: www.fishersci.ca