

Rapid and Robust Protein Chromatography

Chromolith® WP 300 RP-18 2 mm I.D. HPLC Columns

The most hydrophobic of the Chromolith® WP 300 line, the RP-18 column is useful for the resolution of peptides and smaller proteins. One critical quality attribute (CQA) required by regulatory bodies is the peptide map of a biotherapeutic. Peptide maps generated by RP-HPLC provide valuable information about protein structure, stability, and purity. To be effective, the RP-HPLC column must be able to resolve a high percentage of the peptides in the sample. The more peptides, the better the information. The Chromolith® WP 300 RP-18 column gives unsurpassed RP-HPLC resolution of peptide maps from enzymatic digests. The improvements in silica and bonded-phase chemistry incorporated into the Chromolith® WP 300 line improve resolution by increasing efficiency and reducing peak tailing.

Key Benefits:


- Monolithic skeleton permits high flow rates to maximize throughput
- 300 Å mesopores permit large molecules to enter without fear of size-exclusion effects
- Matrix-loaded samples can be injected onto the column with little to no sample prep.

Protein Analysis on Chromolith® WP 300 RP-18

For large molecule separations, high efficiency separations are necessary in order to achieve resolution and good peak shape. Moving to sub-2 µm FPP-packed columns or 2.0 µm SPP-packed columns can deliver that desire; however, this comes at the cost of elevated backpressure. Chromolith® WP 300 RP-18, 2 mm I.D. columns provide UHPLC efficiencies, but at nearly 1/10th the backpressure.

Chromatographic conditions:

Columns:	Chromolith® WP 300 RP-18 100-2 mm (1.52370.0001) SPP, C18, 160 Å, 2.0 µm, 100-2.1 mm FPP, C18, 300 Å, 1.7 µm, 100-2.1 mm
Mobile phase:	A: water (0.1% TFA) B: acetonitrile (0.08% TFA)
Gradient:	4% B to 60% B in 10 minutes
Flow rate:	0.38 mL/min
Detection:	UV, 220 nm
Column temperature:	30 °C
Injection volume:	0.5 µL
Sample:	HPLC Protein Mix 1 mg/mL, water 1) Ribonuclease 2) Cytochrome C 3) Holo-Transferrin 4) Apomyoglobin

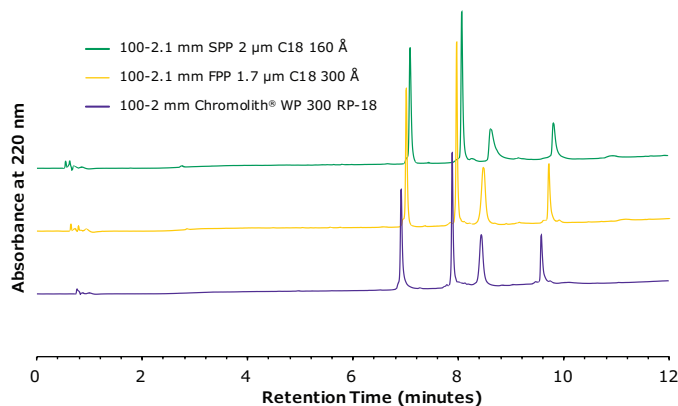


Excellent Lot-to-Lot Reproducibility

Chromolith® WP 300 RP-18 columns exhibit excellent batch-to-batch reproducibility, as demonstrated below with the same peptide map generated for Cytochrome C using three different Chromolith® WP 300 RP-18 columns across three different batches.

Chromatographic conditions:

Column	Chromolith® WP 300 RP-18 100-2 mm (1.52370.0001)		
Mobile Phase	A: acetonitrile 0.08% (v/v) TFA B: water 0.1% (v/v) TFA		
Gradient:	Time	%A	%B
	0	5	95
	25.0	30	70
	30.0	30	70
Flow rate:	0.190 mL/min		
Pressure:	18 bar		
Detection:	Vanquish DAD 20 Hz, UV, 214 nm		
Detector cell:	LightPipe 10 mm		
Temperature:	30 °C		
Injection volume:	0.2 µL		
Sample:	Rapid Trypsin Digestion with SOLu-Trypsin Rapid Digestion Kit 2.5 mg Cytochrome C was added in a PCR vial and dissolved in 320 µL Rapid Trypsin Digestion Buffer. In the solution was 80 µL SOLu-Trypsin added and incubated at 60 °C for 1 hour in a Thermomixer. The digestion was quenched by adding 12 µL of hydrochloric acid 32 %.		

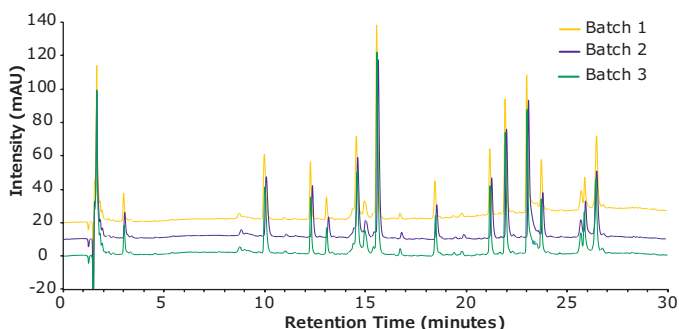


Antibody Fragment Analysis

Fragment analysis of a monoclonal antibody (mAb), also called middle-up analysis, is a useful technique in characterizing mAb domains without the inherent complexity of a peptide map. High efficiency is needed here to resolve subtle, structural variants of the mAb domains. The Chromolith® WP 300 RP-18 column is able to achieve the same separation efficiency and sensitivity as sub-2 µm FPP and 2.0 µm SPP-packed columns but at only 20% of the backpressure of those columns.


Chromatographic conditions:

Column	Chromolith® WP 300 RP-18, 2 mm I.D. (1.52370.0001) SPP, C18, 160 Å, 2.0 µm, 100-2.1 mm FPP, C18, 300 Å, 1.7 µm, 100-2.1 mm	
Mobile Phase	A: Water (0.1% (v/v) TFA) B: Acetonitrile (0.08% (v/v) TFA)	
Gradient:	Time (Min)	%B
	0	20
	1	20
	9	45
Flow rate:	380 µL/min	
Detection:	UV, 220 nm	
Temperature:	80 °C	
Injection volume:	1.0 µL	
Sample:	SigmaMAB, 2 mg/mL (SiLu™ Lite Universal Antibody)	
DTT digest:	60 µL of 40 mM Dithiothreitol (DTT) solution was added in a PCR vial, 40 µL mAb was added and incubated at 37 °C for 30 minutes creating light chain (LC) and heavy chain (HC) parts of the antibody.	

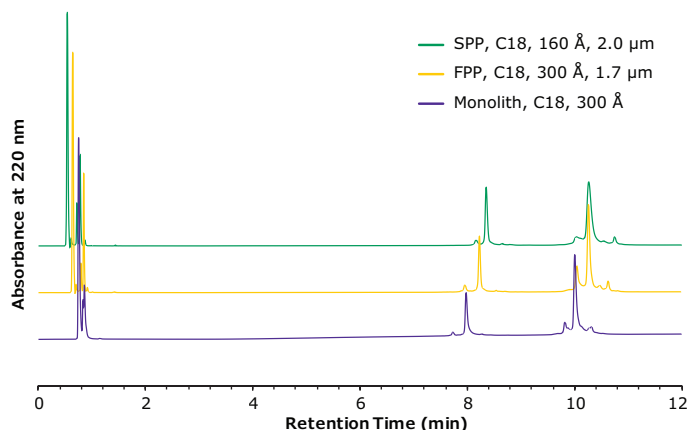

Ordering Information

Cat. No.	Description	Length (mm)	I.D. (mm)
11-010-277	Chromolith® WP 300 RP-18 Column	100	2
11-010-278	Chromolith® WP 300 RP-18 Column	50	2
11-010-279	Chromolith® WP 300 RP-18 Guard Columns (3 units)	5	2

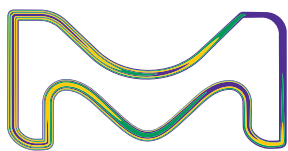

MilliporeSigma is the U.S. and Canada Life Science business of Merck KGaA, Darmstadt, Germany.

© 2022 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved. MilliporeSigma, the vibrant M, Chromolith and Supelco are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.

Lit. No. MS_AD9093EN Ver. 1.0 36577 02/2022

Distributed by Fisher Scientific. Contact us today:

In the United States

Order online: fishersci.com
Call customer service: 1-800-766-7000

In Canada

Order online: fishersci.ca
Call customer service: 1-800-234-7437

